

## Diphenhydramine - pH 9.5, LC-MS

---

Waters Corporation




This is an Application Brief and does not contain a detailed Experimental section.

### Abstract

This application brief demonstrates analysis of diphenhydramine by LC-MS.

## Introduction

The compound analyzed in this study is diphenhydramine.

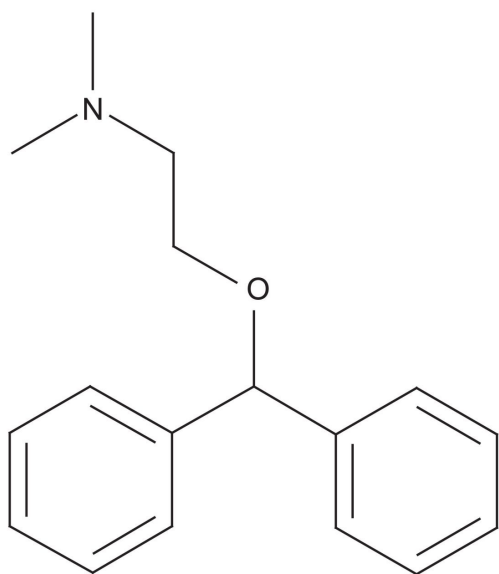

# Diphenhydramine

## Experimental

### Conditions

Column:	Xterra MS C <sub>18</sub> 2.1 x 30 mm, 3.5 μm
Part number:	186000398
Mobile phase A:	0.1% NH <sub>4</sub> OH in H <sub>2</sub> O
Mobile phase B:	0.1% NH <sub>4</sub> OH in ACN

---

Flow rate:	0.2 mL/min to MS
Isocratic mobile phase composition:	55% A; 45% B
Injection volume:	20 $\mu$ L of 100 pg/ $\mu$ L
Temperature:	Ambient
Detection:	MS ESI <sup>+</sup> , SIR 256.12
Instrument:	Alliance 2795 HT, Micromass ZQ

## MS Conditions

Micromass ZQ ESI<sup>+</sup>

Capillary (KV):	3.0
Cone (V):	15
Extractor:	3.0
RF lens:	0.5
Source temp.:	150
Desolvation temp.:	350
Cone gas flow (L/Hr):	60
Desolvation gas flow (L/Hr):	500
LM resolution:	15


Micromass ZQ ESI<sup>+</sup>

HM resolution: 15

Ion energy: 1.0

Multiplier (V): 650

## Results and Discussion

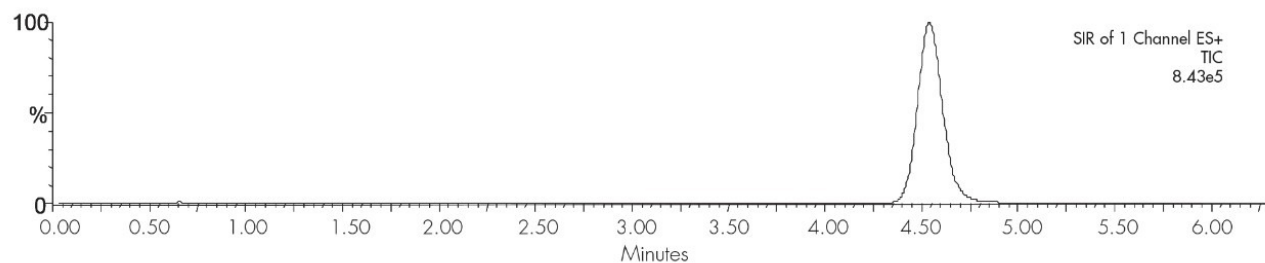

## Featured Products

· [Alliance HPLC System <https://www.waters.com/534293>](https://www.waters.com/534293)

WA20738.040, June 2002



© 2021 Waters Corporation. All Rights Reserved.