


Nota de aplicación

## DNPH Derivatives - pH 2.45

---

Waters Corporation

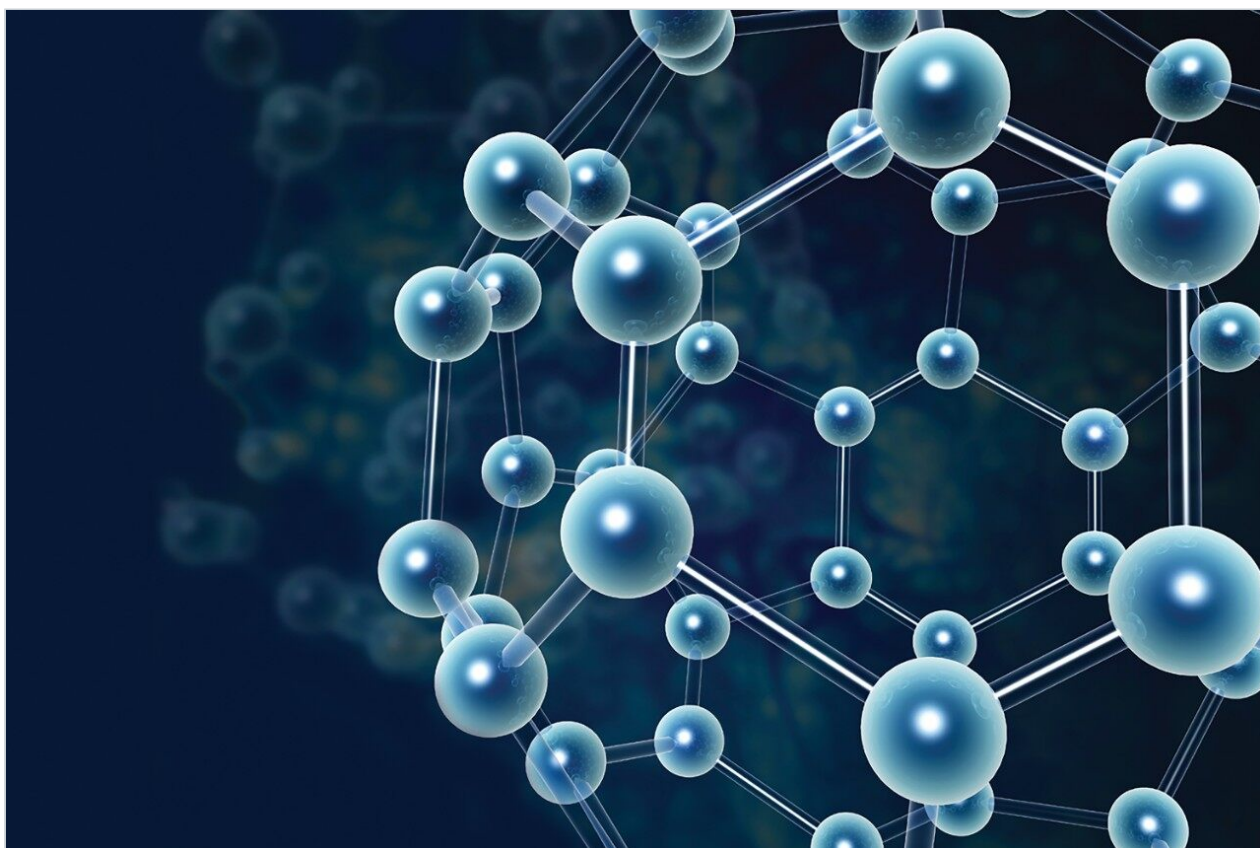

This is an Application Brief and does not contain a detailed Experimental section.

---

### Abstract


This application brief demonstrates analysis of DNPH derivatives.

---


## Introduction

The compounds used in this study are –


1. Formaldehyde-DNPH
2. Acetaldehyde-DNPH
3. Acetone-DNPH
4. Crotonaldehyde-DNPH
5. Cyclohexanone-DNPH

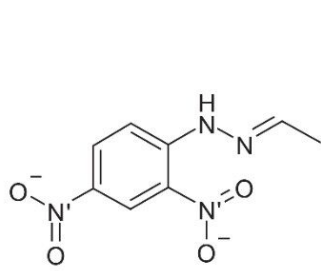


Acetaldehyde-DNPH

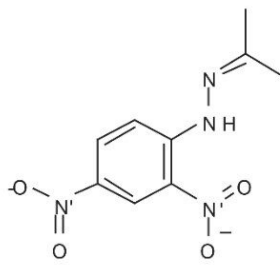


Acetone-DNPH

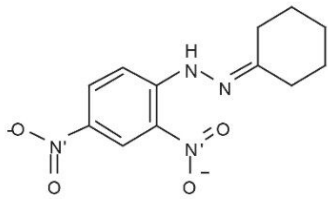

Cyclohexanone-DNPH

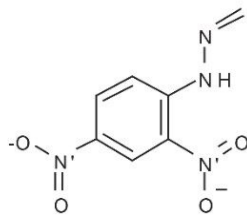

Formaldehyde-DNPH

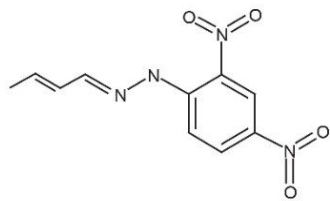

Crotonaldehyde-DNPH

---

## Experimental

### Conditions

Column:

Xterra Phenyl, 4.6 x 150 mm, 5  $\mu$ m

Part number: 186001146

Mobile phase A: H<sub>2</sub>O

Mobile phase B: ACN

Mobile phase C: 50 mM HCOOH, pH 2.45

Flow rate: 1.4 mL/min

Injection volume: 10 µL

Temperature: 30 °C

Detection: UV @ 254 nm


Instrument: Alliance 2695, 2996 PDA

## Gradient

Time (min)	Profile		
	%A	%B	%C
0.0	40	50	10
4.0	40	50	10
10.0	0	90	10

---

## Results and Discussion

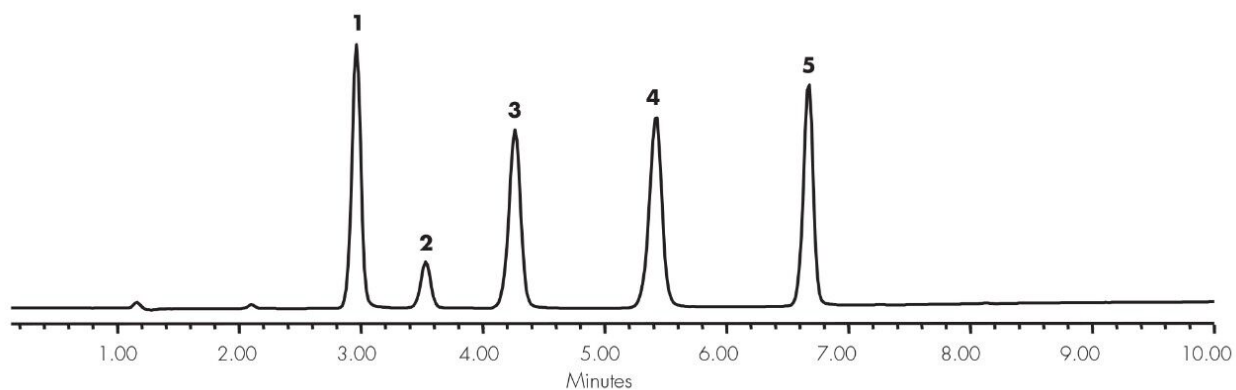

---

## Featured Products

Alliance HPLC System <<https://www.waters.com/534293>>

WA20738.043, June 2002