


RP-HPLC Method for Purification of Dually-Labeled Synthetic Oligonucleotides

Waters Corporation

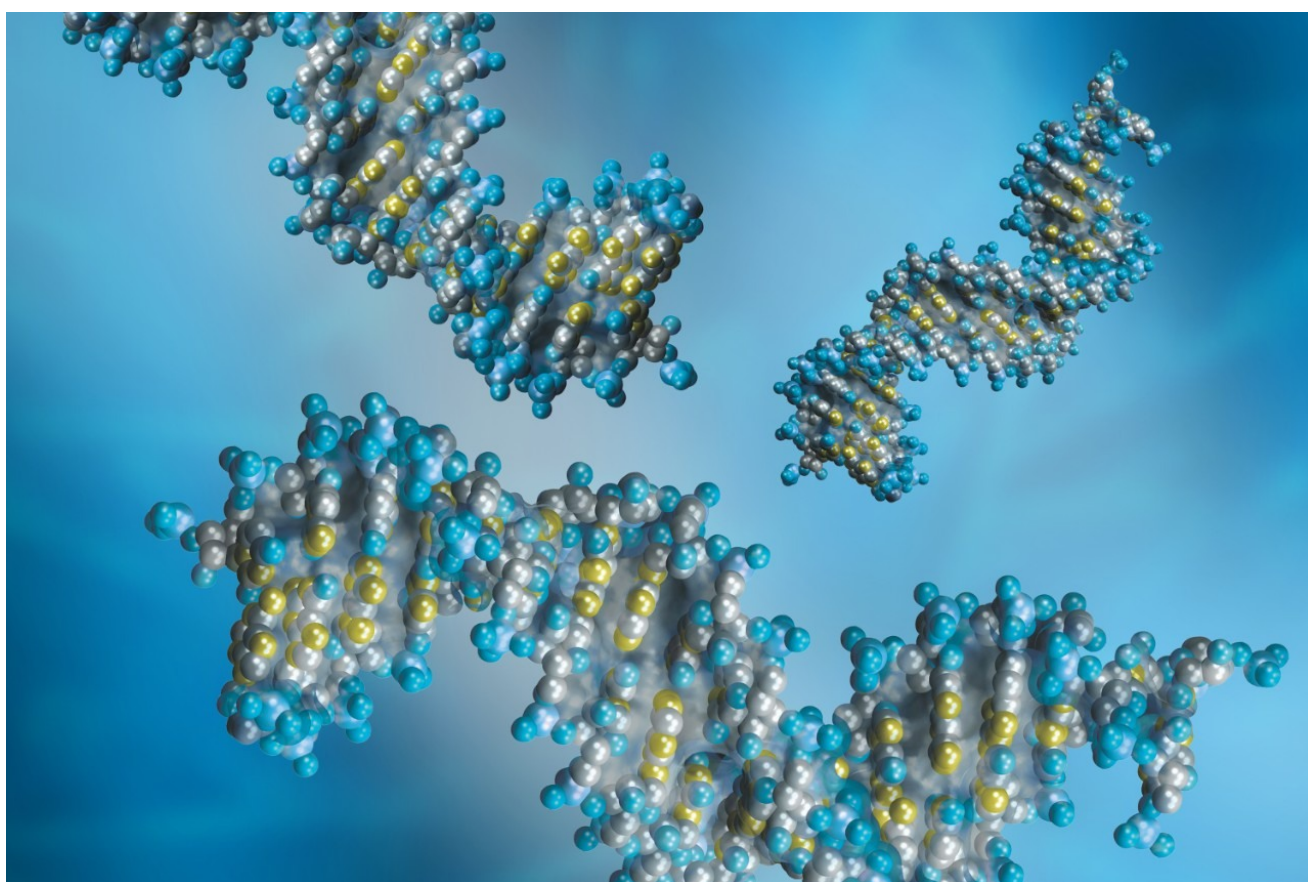

Abstract

This application note demonstrates RP-HPLC method for purification of dually-labeled synthetic oligonucleotides.

Introduction

Dually-Labeled Oligonucleotide Probes

Dually-labeled oligonucleotides are used mainly in quantitative polymerase chain reactions (qPCR), genotyping, and for diagnostic applications. Most common dually-labeled oligonucleotides are TaqMan probes, labeled with 5'FAM and 3' TAMRA dyes. Synthesis of these oligonucleotides is challenging, especially when a "one-pot" approach is adopted without semi-product purification. A target product is contaminated with the partially labeled, non-labeled oligonucleotides, and excess of dyes. The impurities interfere with molecular biology assays and have to be re-moved.

Results and Discussion

XTerra Column for Fast and High-Performance TaqMan Probes Purification

XTerra MS C₁₈ Columns are packed with porous 2.5 µm, hybrid particles. The sorbent has extended stability at temperatures and pH's typically used for oligonucleotide separations (50-60 °C; pH 7-9). Non-labeled oligonucleotides, singly-labeled failure sequences, and excess of free dyes can be easily removed from the target product using XTerra MS C₁₈ Columns. Figure 1 shows the purification of 100 nmole of 31mer TaqMan oligonucleotide produced in a "one-pot" synthesis. Undesirable failed products can be can be successfully removed from product using an XTerra Column. The mobile phases are volatile and can be removed from collected fractions by evaporation. The 4.6 x 50 mm column is suitable for purification of 50–200 nmole scale synthesis in a single injection. Larger amounts (200 nmole–1 µmole) of oligonucleotides can be purified using 10 x 50 mm column.

HPLC system: Alliance® 2695 Separations Module
Column: XTerra®MS C₁₈, 2.5 μm, 4.6 x 50 mm
Part Number: 186000602
Flow rate: 0.5 ml/min
Solvent A: 5 % ACN in 0.1 M TEAA, pH 7
Solvent B: 40 % ACN in 0.1 M TEAA, pH 7

Gradient: From 17.1 % B to 60% B in 30 min
Injection volume: 5-50 μl (5-50 nmole of oligonucleotide)
Column temp.: 60 °C
Detection: UV @ 260 nm and real time UV-VIS 220-600 nm


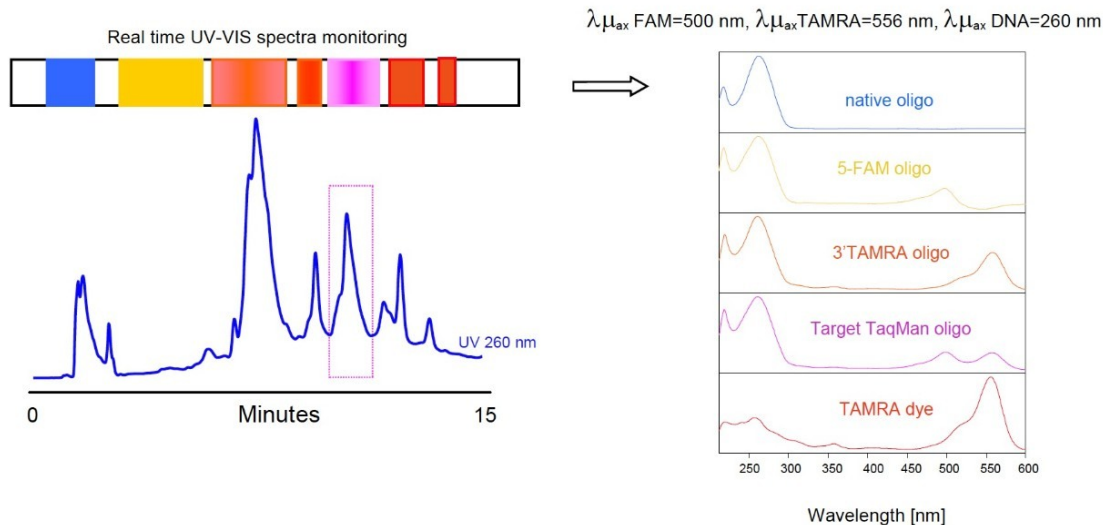


Figure 1. HPLC purification of 100 nmole of 36mer TaqMan oligonucleotide synthesized using “one-pot” approach.

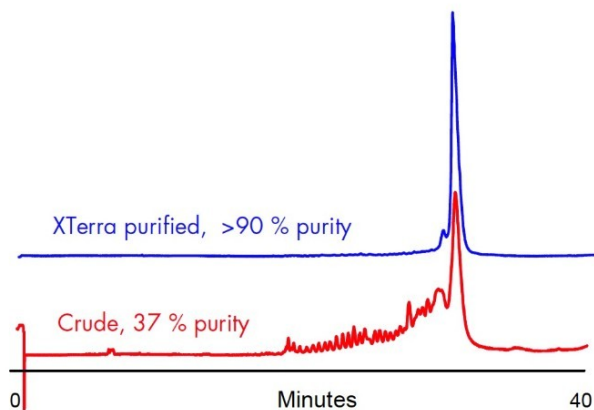

CGE system: Beckman-Coulter P/ACE™ MDQ
Column: BioRad BioCAP™ 75 μm, 34.5 cm
Separation: 15 kV
Sieving matrix: PEG 35 kDa, 2 g/10 ml 0.1 mM Tris-Boric acid, pH 8.3
Injection: 4 s @ 5kV
Column temp.: 30 °C
Detection: UV @ 260 nm

Oligonucleotide was purified on XTerra column (Figure 1)

Figure 2. Capillary Gel Electrophoresis (CGE) quality control of 36mer TaqMan oligonucleotide.

HPLC Purification Strategy

Retention of dye-labeled oligonucleotides is governed by the hydrophobicity the dye. This is demonstrated on the purification of three different TaqMan probes (Figure 3). Non-labeled oligonucleotides elute first (1), followed by FAM labeled failed sequences (2). TAMRA-only labeled failure fragments (3) elute before the dually-labeled target product (4). Other peaks are mostly un-conjugated TAMRA dye. Target product fraction appears to be a group of co-eluting peaks. The partial separation is caused by the differences in HPLC retention of dye

isomers. Mass spectrometry analysis confirmed that the collected product is homogeneous (one molecular mass).

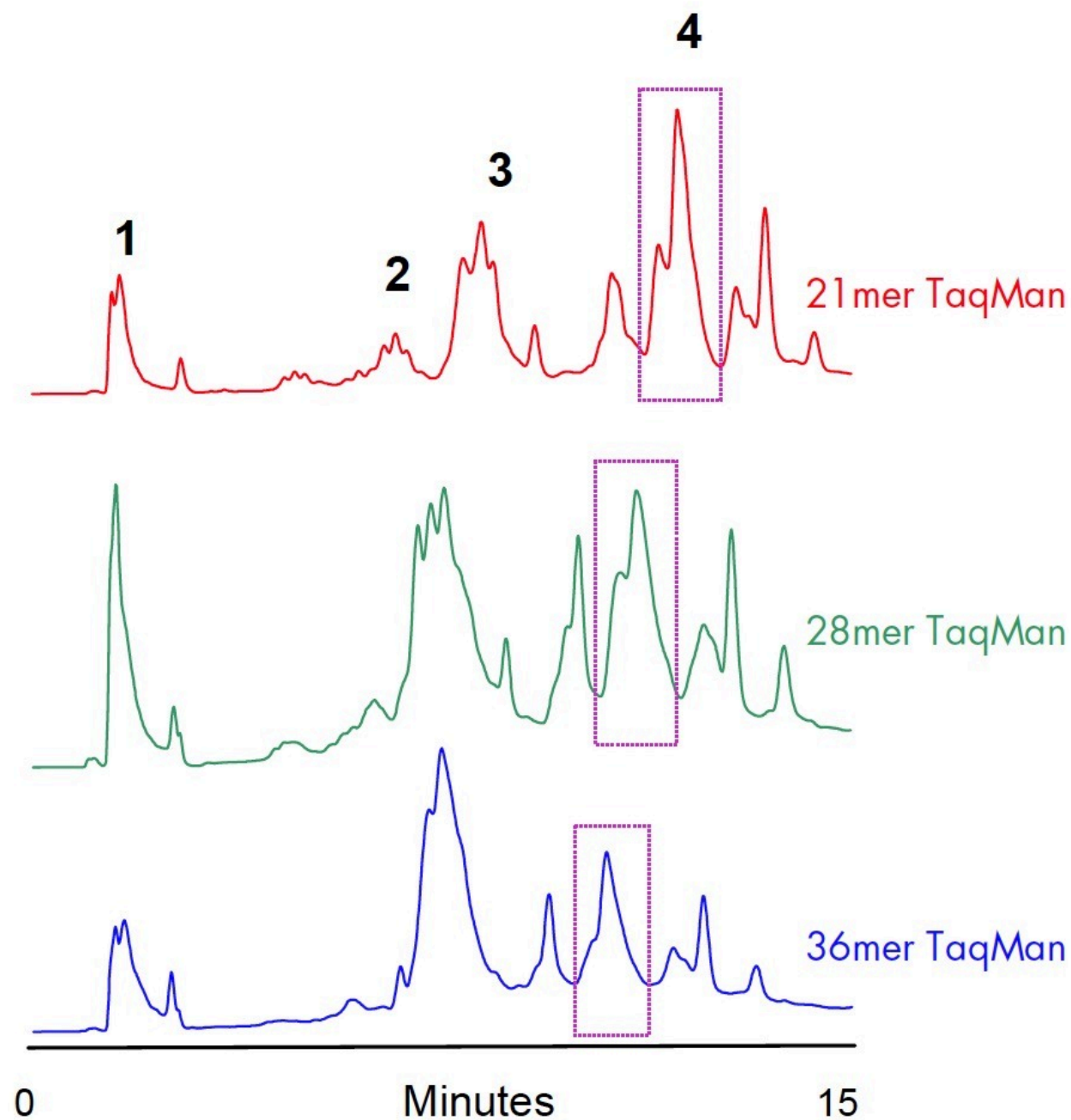


Figure 3. HPLC purification of TaqMan Probes.

In order to select the collection time window, we utilized real-time UV-VIS spectra monitoring along with the 260 nm UV detection (Figure 1). The typical maxima at UV-VIS spectra indicated the elution of non-labeled, singly

labeled and dually-labeled products (three λ_{max} were present in the spectrum). Because of the similar pattern of elution, the purification of different length TaqMan probes becomes a routine for experienced user. Triethylammonium acetate/ACN buffer is volatile. Collected fractions are simply dried down and oligonucleotides are ready for assays.

Featured Products

- [Alliance HPLC System <https://www.waters.com/534293>](https://www.waters.com/534293)

WA20772, June 2003



© 2021 Waters Corporation. All Rights Reserved.