

Chemical Stability Study of ACQUITY UPLC BEH Amide Columns

Waters Corporation


This is an Application Brief and does not contain a detailed Experimental section.

Abstract


This application brief highlights the Chemical stability study of ACQUITY UPLC BEH Amide Columns.

Introduction


Structures

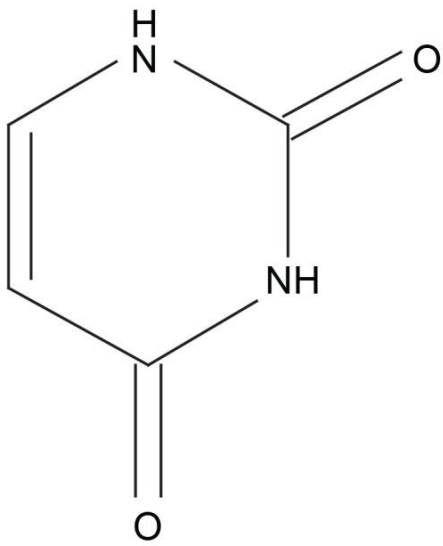

Uracil

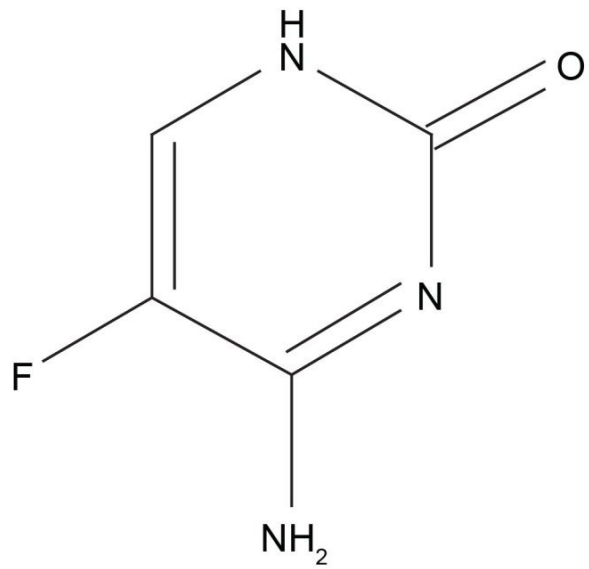

5-Fluorocytosine

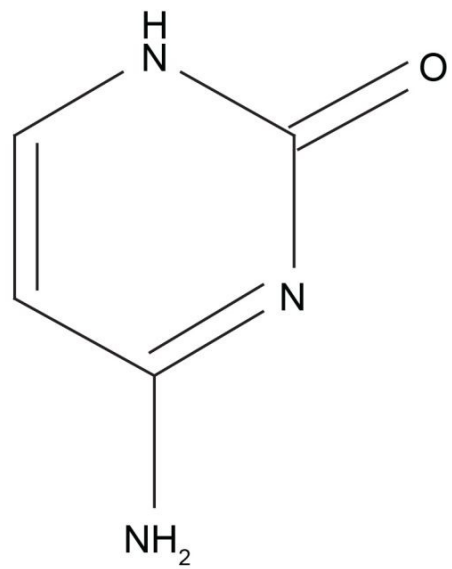

Cytosine

Experimental

Test Conditions

| | |
|-----------------------|--|
| Columns: | ACQUITY UPLC BEH Amide, 2.1 x 50 mm, 1.7 μ m |
| Part Number: | 186004800 |
| Mobile Phase A: | 50/50 MeCN/H ₂ O with 10 mM CH ₃ COONH ₄ , pH 5.5 |
| Mobile Phase B: | 95/5 MeCN/H ₂ O with 10 mM CH ₃ COONH ₄ , pH 5.5 |
| Flow Rate: | 0.5 mL/min |
| Injection Volume: | 2.0 μ L (full loop injection mode) |
| Sample Concentration: | 25 μ g/mL each |
| Sample Diluent: | 75/25 MeCN/MeOH |
| Column Temperature: | 30 °C |
| Weak Needle Wash: | 95/5 MeCN/H ₂ O |
| Detection: | UV @ 254 nm |
| Sampling Rate: | 40 pts/sec |
| Filter Time Constant: | 0.1 |
| Instrument: | Waters ACQUITY UPLC with ACQUITY UPLC PDA Detector |


Gradient

| Time (min) | Profile | |
|------------|---------|----|
| | %A | %B |
| Initial | 1 | 99 |
| 2.00 | 99 | 1 |
| 2.10 | 1 | 99 |
| 2.50 | 1 | 99 |

Results and Discussion

The compounds used in this study are:

1. Uracil
2. 5-fluorocytosine
3. Cytosine

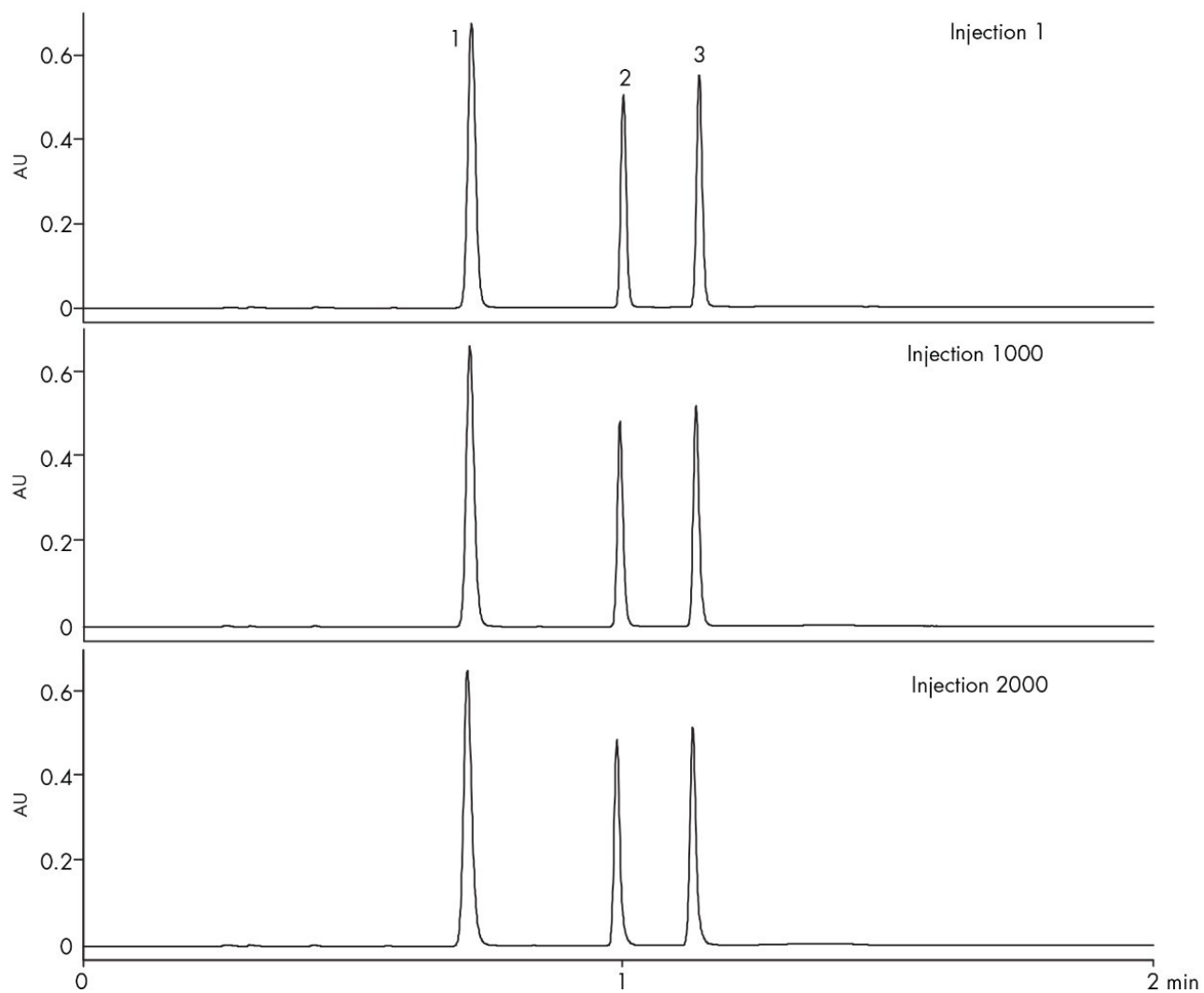

Featured Products

ACQUITY UPLC PDA Detector <<https://www.waters.com/514225>>

WA60106, June 2009

© 2022 Waters Corporation. All Rights Reserved.