

Waters™

應用手冊

Streamlining Analysis of Impurities in the Pharmaceutical Products using Empower 3 ICH Impurity Processing

Margaret Maziarz, Mark Wrona, Paul D.Rainville, Mia Summers, Jade Byrd

Waters Corporation



This is an Application Brief and does not contain a detailed Experimental section.

Abstract

This application brief demonstrates the use of Empower 3 ICH Impurity Processing for streamlining the analysis of impurities of Ziprasidone HCl by defining allowable threshold limits and quickly identifying results above the limits.

Benefits

Utilizing the Empower 3 ICH Impurity Processing enables users to define allowable threshold limits for impurities and quickly identify whether results pass these limits.

Introduction

Impurity profiling including identity and quantity in the drug substances or drug products are a requirement with which every manufacturer must comply. Impurities that develop from the active pharmaceutical ingredient (API) during the formulation and development process of drug product need to be assessed quickly and accurately. Presence of the impurities can compromise safety and efficacy of the end pharmaceutical product and must be effectively monitored. The International Committee for Harmonization (ICH) has published guidelines on impurities in drug substances¹ and drug products,² providing allowable threshold limits for impurities to monitor safety.

The Empower 3 ICH Impurity Processing function simplifies quantitative analysis by quickly identifying impurities above the ICH allowable limits defined by the user. Empower Software automatically compares the calculated amount of impurities against the limits and flags any failing results. This allows quick evaluation of the pharmaceutical products for safety during formulation or release testing.

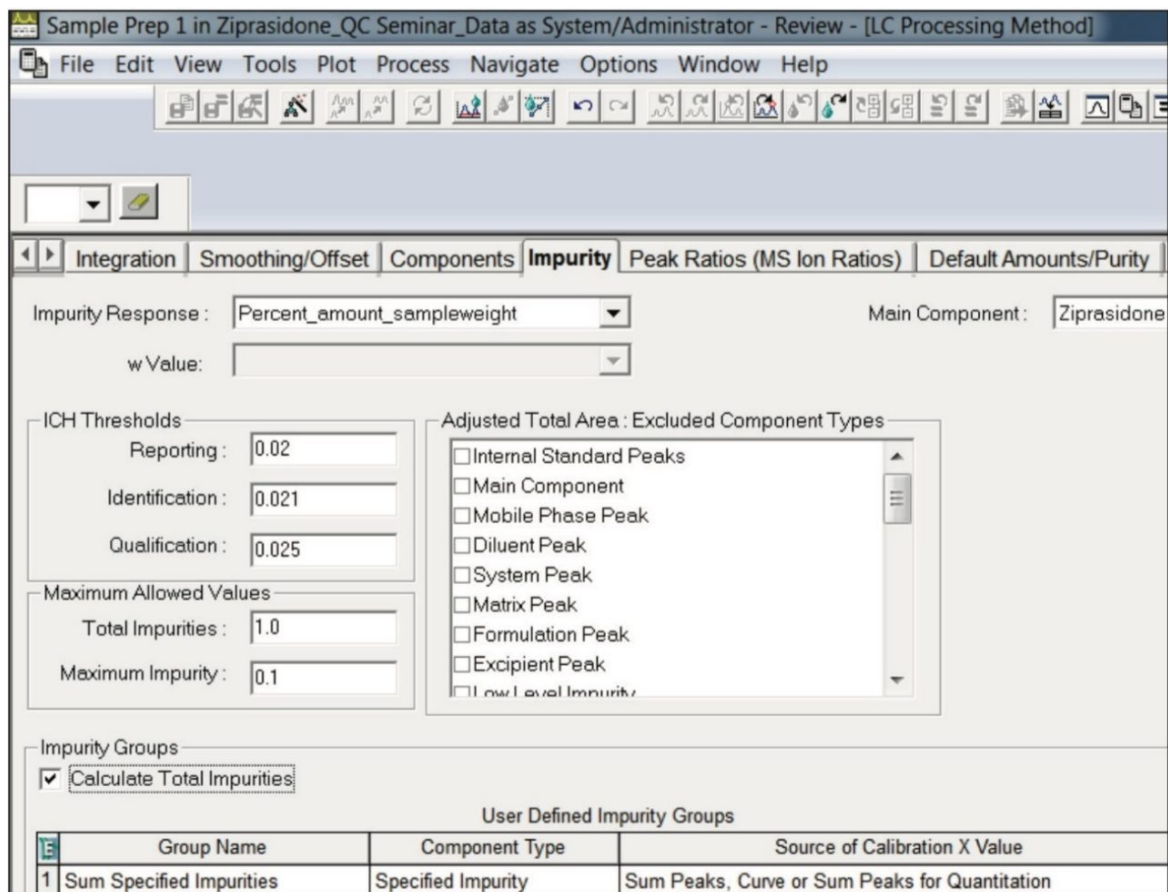


Figure 1. Empower 3 ICH Impurity Processing. Selection of field for impurity response, ICH thresholds, maximum allowed values, and group name to sum specified impurities.


This technology brief illustrates use of Empower 3 ICH Impurity Processing for streamlining quantitative analysis of impurities of Ziprasidone HCl in a capsule formulation.

Results and Discussion

The UPLC method for Ziprasidone HCl and its USP-specified related substances was previously developed in the Waters application note 720004639EN.³ Empower 3 ICH Impurity Processing was used to streamline the quantitative analysis of related substances in the Ziprasidone HCl capsule formulation by defining the ICH threshold limits and identifying results that did not pass these limits. First, we specified the formula for calculating quantity of impurities or impurity response in the Empower processing method (Figure 1). Next, we defined the ICH thresholds for reporting, identification, and qualification of impurities. Then, we specified maximum allowed values for the total impurities and maximum impurity. In this example, we selected more

stringent criteria compared to the ICH guidelines to demonstrate the workflow of identifying and flagging peaks that exceed impurity threshold limits. We used a reporting threshold of 0.02% of API. This is lower than the ICH guidelines for reporting impurities in new drug products of 0.1% based on maximum daily dosage of ≤ 1 g.² In addition, we can select tighter limits than the generic ICH thresholds for each specified impurity of the pharmaceutical product. Furthermore, we can group specific types of impurities to calculate their total.

Finally, we processed the chromatographic data to determine quantity of impurities. The data (Figure 2A) shows that two impurities, compounds A and B, were detected in the capsule sample, using Empower's proprietary peak detection algorithm ApexTrack. Empower Software quickly identified the out-of-specification (OOS) result for compound A (flagged in red), as this value was above the threshold limits defined in the processing method (Figure 2B). The results for analysis of impurities can be easily reported (Figure 3). The report template is customizable to tailor the reporting requirements for each user's needs and can facilitate the use of electronic signatures as shown in Figure 3.

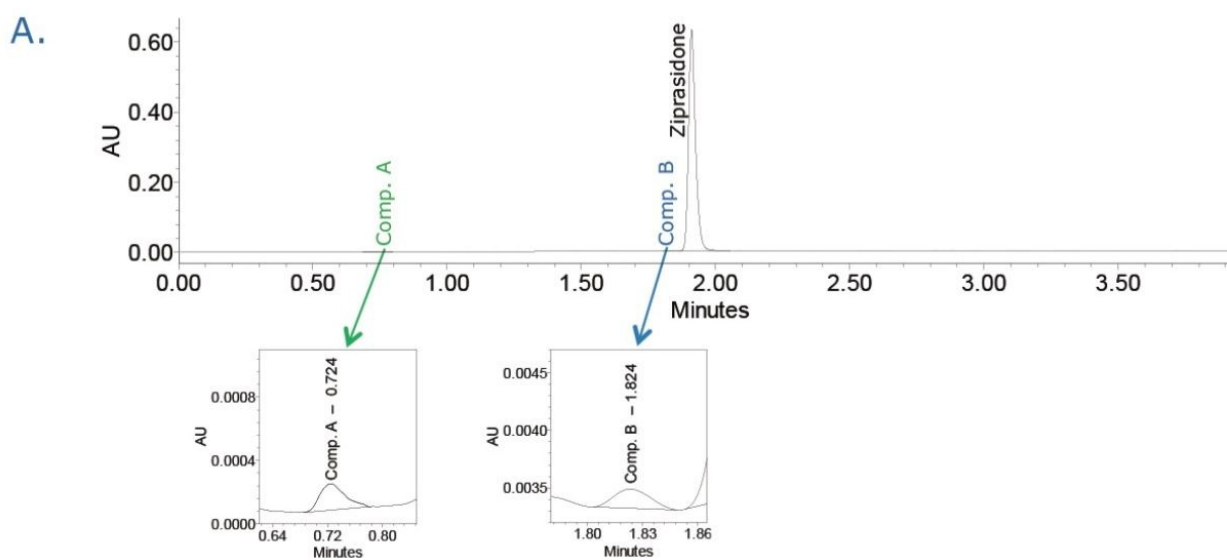

B.

Name	Component Type	Retention Time (min)	Area ($\mu\text{V}\cdot\text{sec}$)	Impurity Response	Above Reporting Threshold	Above Identification Threshold	Above Qualification Threshold	ICH Threshold	Pass?
Compound A	Specified Impurity	0.724	439	0.027	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	<input checked="" type="checkbox"/>	Above Qualification Threshold	Fail
Compound B	Specified Impurity	1.824	230	0.017	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	Below Reporting Threshold	Pass
Ziprasidone	Main Component	1.911	1145524						
Compound D	Low Level Impurity	2.978							
Compound C	Low Level Impurity	3.119							
Sum Specified Impurities	Specified Impurity		669	0.044					Fail

Figure 2. Analysis of impurities of Ziprasidone HCl in the capsule sample: A. Capsule sample with UV at 229 nm., B. Results for quantitative determination of impurities.

SAMPLE INFORMATION

Sample Name: Sample Prep 1	Sample Set Name: Impurities_UPLC BEH
Sample Type: Unknown	Acq. Method Set: UPLC_BEH
Vial: 1:A,6	Processing Method: Impurities_ICH Imp Process_2
Injection #: 1	Channel Name: 229nm
Injection Volume: 1.40 ul	Proc. Chnl. Descr.: ACQUITY TUV ChA
Run Time: 4.0 Minutes	Calibration Id 5954
Result Set Id 5934	

% Impurity

	SampleName	Name	RT (min)	Sample Weight (mg)	Dilution (mL)	Amount (mg)	Reporting Threshold	% Impurity	Pass?
1	Sample Prep 1	Compound A	0.724	1.301	10	0.000348	0.02	0.027	Fail
2	Sample Prep 1	Compound B	1.824	1.301	10	0.000221	0.02	0.017	Pass
3	Sample Prep 1	Ziprasidone	1.911	1.301	10				
Mean									

ICH Thresholds:

Reporting Threshold: 0.02
 Identification Threshold: 0.021
 Qualification Threshold: 0.025

Sign Off Full Name: Jane Smith (Lab_Manager)
 Sign Off Reason: Sign Off Level 1, Reason: Approval

Figure 3. Empower report for analysis of impurities of Ziprasidone HCl in the capsule sample. Results approved and signed off by lab manager via electronic signature (OOS flags shown in red).

Conclusion

Utilizing the Empower 3 ICH Impurity Processing enables users to define allowable threshold limits for impurities and quickly identify whether results are within these limits. The threshold limits defined by the user are clearly displayed by Empower during data review or in a report. This reduces the time and potential errors associated with manual verification and enhances the confidence that any results above the threshold limits are quickly identified.

Overall, Empower 3 ICH Impurity Processing is a powerful (yet simple) tool that can be utilized by any laboratory to monitor levels of impurities during formulation or release testing of the pharmaceutical products and streamline their quantitative analysis protocols.

References

1. ICH Q3A(R2), Impurities in New Drug Substances, International Conference on Harmonization, October 2006.
2. ICH Q3B(R2), Impurities in New Drug Products, International Conference on Harmonization, June 2006.
3. Maziarz M, Jones MD, Potts WB, Forini F. Transfer of Two UPS Compendial Methods for Impurities of Ziprasidone HCl to a Single UPLC Method. Waters Application Note 720004639EN, April 2013.

Featured Products

- [Empower 3 Chromatography Data Software <https://www.waters.com/10190669>](https://www.waters.com/10190669)

720005180, September 2014



©2019 Waters Corporation. All Rights Reserved.