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
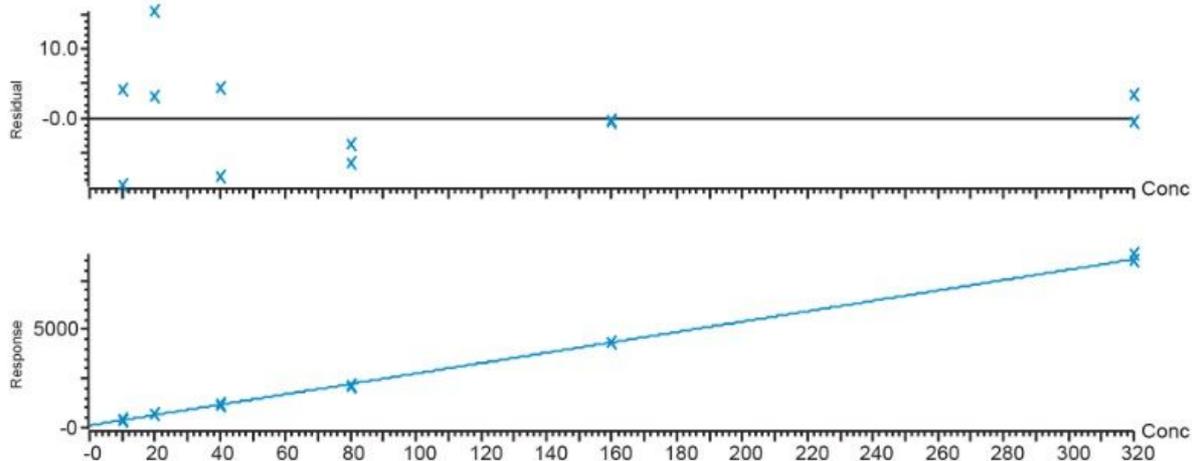
Compound name: 17 Alpha Ethynyl Estradiol  
 Correlation coefficient:  $r = 0.999001$ ,  $r^2 = 0.998004$   
 Calibration curve:  $26.4155 * x + 120.806$   
 Response type: External Std, Area  
 Curve type: Linear, Origin: Exclude, Weighting: 1/x, Axis trans: None

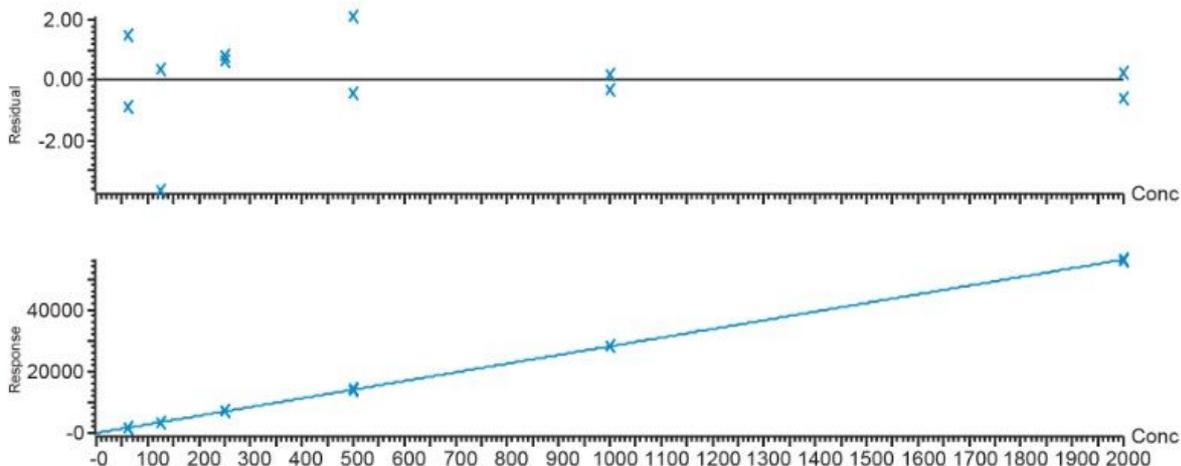Compound name: 17 Beta Estradiol  
 Correlation coefficient:  $r = 0.999950$ ,  $r^2 = 0.999901$   
 Calibration curve:  $28.1779 * x + -59.3078$   
 Response type: External Std, Area  
 Curve type: Linear, Origin: Exclude, Weighting: 1/x, Axis trans: None

Compound name: Estrone  
 Correlation coefficient:  $r = 0.999910$ ,  $r^2 = 0.999819$   
 Calibration curve:  $171.91 * x + 336.28$   
 Response type: External Std, Area  
 Curve type: Linear, Origin: Exclude, Weighting: 1/x, Axis trans: None

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Compound name: 17 Alpha Ethynyl Estradiol  
Correlation coefficient:  $r = 0.998732$ ,  $r^2 = 0.997467$   
Calibration curve:  $17.5798 * x + 296.967$   
Response type: External Std, Area  
Curve type: Linear, Origin: Exclude, Weighting:  $1/x$ , Axis trans: None  
Standard Addition Concentration : 16.8925

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---

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1. Schwindt A, Winkelman D, Keteles L, *et al.* An Environmental Oestrogen Disrupts Fish Population Dynamics through Direct and Transgenerational Effects on Survival and Fecundity. *Journal of Applied Ecology*. 2014.
2. Directive 2013/39/EU of the European Parliament and of the Council of 12 August 2013 Amending Directive 2000/60/EC and 2008/105/EC in Respect of Priority Substances in the Field of Water Policy Text with EEA Relevance. August 2013.
3. Directive 2000/60/EC of the European Parliament and of the Council of 23 October 2000 Establishing a Framework for Community Action in the Field of Water Policy. October 2000.
4. Commission Implementing Decision (EU) 2015/495 of 20 March 2015 Establishing a Watch List of Substances for Union-Wide Monitoring in the Field of Water Policy Pursuant to Directive 2008/105/EC of the European Parliament and of the Council (Notified under Document C(2015) 1756) Text with EEA Relevance. March 2015.
5. Ross E, Boag A, Todd H, Gatward N. Analysis of Natural and Synthetic Estrogens at Sub-PPT Levels in Surface Water and Crude Influent Water Utilizing the ACQUITY UPLC System with 2D LC Technology and Xevo TQ-S. Waters Technology Brief No. 720005626en <<https://www.waters.com/nextgen/us/en/library/application-notes/2016/analysis-of-natural-and-synthetic-estrogens-at-sub-ppt-in-water-lcmsms.html>> . March 2016.
6. Chang H. *et al.* Occurrence of Androgens and Progestogens in Wastewater Treatment Plants and Receiving River Waters: Comparison to Estrogens. *Waters Research*. 45(2):732–740. 2011.

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TargetLynx <<https://www.waters.com/513791>>

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©2019 Waters Corporation. All Rights Reserved.