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1. Baksam V, Nimmakayala S, Devineni SR, Muchumarri RM, Shandilya S, Kumar P. Isolation and Characterization of Thermal Degradation Impurity in Brimonidine Tartrate by HPLC, LC-MS/MS, and 2DNMR. *Journal of Pharmaceutical and Biomedical Analysis*.2021.(205) 114297.
2. Aliyu AO, Garba S, Balogun LO, Awe FE. Quality Assessment of Some Selected Brands of Amoxicillin Clavulanate from Pharmaceutical Stores in Kaduna Metropolis, Nigeria. *Journal of Pharmacy and Biological Sciences*.2021-Apr.(16) 33-40.
3. Monisha ST, Ela KN, Islam R, Ether SA, Rahman FI. Quality Attributes Comparison of Selected Brands of Rosuvastatin Calcium Tablets Marketed in the US and Bangladesh. *Journal of Pharmaceutical Research International*. 2021 (33) 46-55.
4. 621-Chromatography.pdf Accessed 9-Sept-2021. <https://www.bioglobax.com/wp-content/uploads/2018/08/621-Chromatography.pdf> <<https://www.bioglobax.com/wp-content/uploads/2018/08/621-Chromatography.pdf>>.

5. Swann T, Nguyen JM. USP Method Modernization Using “Equivalent L/dp” and “Equivalent N” Allowed Changes with CORTECS C₈ and CORTECS UPLC C₈ Columns. Waters Application Note. 2016, [720005666EN](#).
6. Naproxen-sodium-tablets.pdf. Naproxen Sodium Tablets USP monograph. Accessed 9-Sept-2021.
https://www.uspnf.com/sites/default/files/usp_pdf/EN/USPNF/iras/naproxen-sodium-tablets.pdf <
https://www.uspnf.com/sites/default/files/usp_pdf/EN/USPNF/iras/naproxen-sodium-tablets.pdf>.
7. Berthelette KD, Summers M, Fountain KJ. Improving Resolution using CORTECS UPLC Columns. Waters Application Note. 2013, [720004737EN](#).
8. Berthelette KD, Nguyen JM, Turner JE. Improving Peak Capacity while Maintaining Selectivity using CORTECS Columns on an Agilent 1290 LC system. Waters Application Note. 2021, [720007250EN](#).
9. Shia JC, Yabu M, Tran KV, Young M. Improved Resolution of Paraquat and Diquat: Drinking Water Analysis Using the CORTECS UHPLC HILIC Column. Waters Application Note. 2013, [720004732EN](#).
10. Danaceau JP, Chambers EC, Fountain KJ. Advantages of CORTECS C₁₈ 2.7 μm and XBridge BEH Phenyl XP 2.5 μm Columns for the Analysis of a Comprehensive Panel of Pain Management Drugs for Forensic Toxicology. Waters Application Note. 2014, [720005185EN](#).
11. Berthelette KD, Fountain KJ, Summers M. Improving Method Robustness for Routine Analysis of Pharmaceutical Formulations. Waters Application Note. 2013, [720004807EN](#).
12. Acetonitrile Suitable for HPLC, gradient grade ≥99.9% Product page. Cost of 1L: \$185.00.
<https://www.sigmaaldrich.com/US/en/product/sigald/34851?context=product> <
<https://www.sigmaaldrich.com/US/en/product/sigald/34851?context=product>>.
13. Water Suitable for HPLC Product page. Cost of 1 L: \$51.20.
<https://www.sigmaaldrich.com/US/en/product/sigald/270733?context=product> <
<https://www.sigmaaldrich.com/US/en/product/sigald/270733?context=product>>.
14. Acetic Acid, Glacial HPLC, Meets ACS Specifications Product Page. Cost of 500 mL: \$87.90.
<https://www.sigmaaldrich.com/US/en/product/mm/ax0074?context=product> <
<https://www.sigmaaldrich.com/US/en/product/mm/ax0074?context=product>>.
15. Calculation of cost per L of mobile phase.

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© 2022 Waters Corporation. All Rights Reserved.