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1. Centers for Disease Control and Prevention.What is Diabetes?. Centers for Disease Control and Prevention.2021 [cited 2022 Apr 28].Available from: <https://www.cdc.gov/diabetes/basics/diabetes.html>.
2. Metformin.The Johns Hopkins Patient Guide to Diabetes.*John Hopkins Medicine*; 2016 [cited 2022 Apr 28].Available from: <https://hopkinsdiabetesinfo.org/medications-for-type-2-diabetes-metformin/>.
3. Walter T, Berthelette K, Patel A, Alden B, McLaughlin J, Field J, Lawrence N, Shiner S.Introducing Atlantis BEH Z-HILIC: A Zwitterionic Stationary Phase Based on Hybrid Organic/Inorganic Particles, Waters Application Note, 2021, [720007311](#).
4. Clements B, Maziarz M, Rainville P. Method for the Analysis and Quantitation of Pharmaceutical Counterions Utilizing Hydrophilic Interaction Liquid Chromatography.Waters Application Note, 2022, [720007564](#).
5. Bethesda (MD): National Library of Medicine (US), National Center for Biotechnology Information; *PubChem*. PubChem Compound Summary for CID 10005, Cyanoguanidine; [cited 2022 Apr. 25].2004. Available from: <https://pubchem.ncbi.nlm.nih.gov/compound/cyanoguanidine#:~:text=Molecular%20Weight,-84.08,-Dates> < <https://pubchem.ncbi.nlm.nih.gov/compound/cyanoguanidine#:~:text=Molecular%20Weight,-84.08,-Dates>> .
6. Bethesda (MD): National Library of Medicine (US), National Center for Biotechnology Information; *PubChem* .PubChem Compound Summary for CID 24196656; [cited 2022 Apr. 25].2004. Available from: <https://pubchem.ncbi.nlm.nih.gov/compound/24196656#:~:text=Reference-,Molecular%20Weight,204.62,-Computed%20by%20PubChem> < <https://pubchem.ncbi.nlm.nih.gov/compound/24196656#:~:text=Reference-,Molecular%20Weight,204.62,-Computed%20by%20PubChem>> .
7. Bethesda (MD): National Library of Medicine (US), National Center for Biotechnology Information; *PubChem* .PubChem Compound Summary for CID 11469186, 1,3,5-Triazine-2,4,6-triamine, N,N'-dimethyl-; [cited 2022 Apr. 25].2004. Available from: https://pubchem.ncbi.nlm.nih.gov/compound/2-N_4-N-dimethyl-1_3_5-triazine-2_4_6-triamine#:~:text=Molecular%20Weight,-154.17,-Computed%20by%20PubChem < https://pubchem.ncbi.nlm.nih.gov/compound/2-N_4-N-dimethyl-1_3_5-triazine-2_4_6-triamine#:~:text=Molecular%20Weight,-154.17,-Computed%20by%20PubChem> .
8. Bethesda (MD): National Library of Medicine (US), National Center for Biotechnology Information; *PubChem* .PubChem Compound Summary for CID 7955, Melamine; [cited 2022 Apr. 25].2004. Available from:

<https://pubchem.ncbi.nlm.nih.gov/compound/melamine#:~:text=Reference-,Molecular%20Weight,126.12,-Computed%20by%20PubChem> <
<https://pubchem.ncbi.nlm.nih.gov/compound/melamine#:~:text=Reference-,Molecular%20Weight,126.12,-Computed%20by%20PubChem>> .

9. Bethesda (MD): National Library of Medicine (US), National Center for Biotechnology Information; *PubChem*. PubChem Compound Summary for CID 15486, 1-Methylbiguanide; [cited 2022 Apr. 25].2004. Available from: <https://pubchem.ncbi.nlm.nih.gov/compound/1-Methylbiguanide#:~:text=More...-,Molecular%20Weight,115.14,-Dates> <<https://pubchem.ncbi.nlm.nih.gov/compound/1-Methylbiguanide#:~:text=More...-,Molecular%20Weight,115.14,-Dates>> .
10. Bethesda (MD): National Library of Medicine (US), National Center for Biotechnology Information; *PubChem*. PubChem Compound Summary for CID 10473, Dimethylamine hydrochloride; [cited 2022 Apr. 25].2004. Available from: <https://pubchem.ncbi.nlm.nih.gov/compound/Dimethylamine-hydrochloride#:~:text=4.2Mass%20Spectrometry> <
<https://pubchem.ncbi.nlm.nih.gov/compound/Dimethylamine-hydrochloride#:~:text=4.2Mass%20Spectrometry>> .

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© 2022 Waters Corporation. All Rights Reserved.

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