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1. UNAIDS. Global HIV & AIDS Statistics — 2020 Fact Sheet. 2021. Available from: <https://www.unaids.org/en/resources/fact-sheet> <<https://www.unaids.org/en/resources/fact-sheet>>
2. Centers for Disease Control and Prevention. HIV Basics. 2019. Available from: <https://www.cdc.gov/hiv/basics/index.html> <<https://www.cdc.gov/hiv/basics/index.html>>
3. Centers for Disease Control and Prevention. AIDS and Opportunistic Infections. 2019. Available from: <https://www.cdc.gov/hiv/basics/livingwithhiv/opportunisticinfections.html> <<https://www.cdc.gov/hiv/basics/livingwithhiv/opportunisticinfections.html>>
4. Pan American Health Organization. Antiretroviral Therapy - PAHO/WHO | 2023. Available from: [https://www.paho.org/en/topics/antiretroviral-therapy#:~:text=Antiretroviral%20therapy%20\(ART\)%20is%20treatment](https://www.paho.org/en/topics/antiretroviral-therapy#:~:text=Antiretroviral%20therapy%20(ART)%20is%20treatment) <[https://www.paho.org/en/topics/antiretroviral-therapy#:~:text=Antiretroviral%20therapy%20\(ART\)%20is%20treatment](https://www.paho.org/en/topics/antiretroviral-therapy#:~:text=Antiretroviral%20therapy%20(ART)%20is%20treatment)>
5. National Institute of Health. Appendix B: Drug Characteristics Tables | NIH. 2023 [cited 2023 Aug 11]. Available from: <https://clinicalinfo.hiv.gov/en/guidelines/hiv-clinical-guidelines-adult-and-adolescent-arv/drug-characteristics-tables?view=full#table7> <<https://clinicalinfo.hiv.gov/en/guidelines/hiv-clinical-guidelines-adult-and-adolescent-arv/drug-characteristics-tables?view=full#table7>>
6. National Institutes of Health. Enfuvirtide - Patient | NIH. 2023. Available from: <https://clinicalinfo.hiv.gov/en/drugs/enfuvirtide/patient> <<https://clinicalinfo.hiv.gov/en/drugs/enfuvirtide/patient>>
7. PubChem. Enfuvirtide. National Center for Biotechnology Information. 2023. Available from: <https://pubchem.ncbi.nlm.nih.gov/compound/Enfuvirtide> <<https://pubchem.ncbi.nlm.nih.gov/compound/Enfuvirtide>>
8. Charbe N, Baldelli S, Cozzi V, Castoldi S, Cattaneo D, Clementi E. Development of an HPLC–UV Assay Method for the Simultaneous Quantification of 9 Antiretroviral Agents in the Plasma of HIV-Infected Patients. *Journal of Pharmaceutical Analysis*. December 2016. 6(6):396–403.
9. D' Avolio A, Siccardi M, Sciandra M, Lorena B, Bonora S, Trentini L, *et al.* HPLC–MS Method for the Simultaneous Quantification of the New HIV Protease Inhibitor Darunavir, and 11 other Antiretroviral Agents in Plasma of HIV-Infected Patients. *Journal of Chromatography B*. November 2007. 859(2):234–40.

10. Birdsall R, Kellet J, Ippoliti S, Qing Yu Y. Increasing Recovery and Chromatographic Performance of “Acidic” Peptides Using Waters ACQUITY Premier Solution. Waters Application Note. [720007173](#). 2021.
11. Buck MA, Olah TA, Weitzmann CJ, Cooperman BS. Protein Estimation by the Product of Integrated Peak Area and Flow Rate. *Analytical Biochemistry*. Nov 1989. [cited 2023 Aug 11]. 182(2):295–9. Available from: <https://www.sciencedirect.com/science/article/abs/pii/0003269789905976?via%3Dihub> <
<https://www.sciencedirect.com/science/article/abs/pii/0003269789905976?via%3Dihub>>
12. U.S. Food and Drug Administration. Reviewer Guidance, Validation of Chromatographic Methods. *FDA U.S. Food and Drug Administration*. 2018. Available from: <https://www.fda.gov/regulatory-information/search-fda-guidance-documents/reviewer-guidance-validation-chromatographic-methods> <
<https://www.fda.gov/regulatory-information/search-fda-guidance-documents/reviewer-guidance-validation-chromatographic-methods>>
13. Instrument Specifications ACQUITY QDa Detector. Waters Corporation; 2016 [cited 2023 Aug 11]. Available from: <https://www.waters.com/webassets/cms/library/docs/720004799en.pdf> <
<https://www.waters.com/webassets/cms/library/docs/720004799en.pdf>>
14. Chang D, Kolis SJ, Linderholm K, Julian TF, R. Nachi, A.M. Dzerk, et al. Bioanalytical Method Development and Validation for a Large Peptide HIV Fusion Inhibitor (Enfuvirtide, T-20) and Its Metabolite in Human Plasma Using LC–MS/MS. *Journal of Pharmaceutical and Biomedical Analysis*. 2005 Jul 1;38(3):487–96.
15. International Conference on Harmonisation of Technical Requirements for Registration of Pharmaceuticals for Human Use Ich Harmonised Tripartite Guideline Validation of Analytical Procedures: Text and Methodology Q2(R1). *ICH*. November 2005 . Available from: <https://database.ich.org/sites/default/files/Q2%28R1%29%20Guideline.pdf> <
<https://database.ich.org/sites/default/files/Q2%28R1%29%20Guideline.pdf>>

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© 2024 Waters Corporation. All Rights Reserved.

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