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1. Schymanski E, Zhang J, Thiessen P, Chirsir P, Kondic, T, Bolton, E. Per- and Polyfluoroalkyl substances (PFAS) in PubChem: 7 Million and Growing. *Environ.Sci.Technol.* 2023 57 (44): 16918–16928
2. US Environmental Protection Agency. *CompTox Chemicals Dashboard*. PFAS Structures in DSSTox (update August 2022). <https://comptox.epa.gov/dashboard/chemical-lists/PFASSTRUCTV5> <
<https://comptox.epa.gov/dashboard/chemical-lists/PFASSTRUCTV5>>
3. US Environmental Protection Agency. Analysis of Per- and Polyfluoroalkyl Substances (PFAS) in Aqueous, Solid, Biosolids, and Tissue Samples by LC-MS/MS. January 2024
4. ASTM D8421–22, Standard Test Method for Determination of Per- and Polyfluoroalkyl Substances (PFAS) in Aqueous Matrices by Co-solvation followed by Liquid Chromatography Tandem Mass Spectrometry (LC/MS/MS), ASTM International, West Conshohocken, PA, 2022, www.astm.org <<http://www.astm.org/>>
5. Liu Y, D’ Agostino LA, Qu G, Jiang G, Martin JW. High-resolution mass spectrometry (HRMS) methods for non-target discovery and characterization of poly- and perfluoroalkyl substances (PFASs) in environmental and human samples. *Trends Anal Chem* 2019, 115420
6. McCullagh M, Hodgkinson M, Dilllon L, van Bevel B, Burgess JA, Jogsten IE. Determination and Characterization of Perfluoroalkyl and Polyfluoroalkyl Substances (PFAS’ s) in Environmental Samples Using UPLC Ion Mobility MS. Waters Application Note [720005090](#). 2014 June.
7. Dodds JN, Hopkins ZR, Knappe DRU, Baker ES. Rapid Characterization of Per- and Polyfluoroalkyl substances (PFAS) by Ion Mobility Spectrometry-Mass Spectrometry (IMS-MS). *Anal Chem* 2020, 92, 6, 4427–4435
8. Organtini KL, Rosnack KJ, Plummer C, Hancock P, Burt O. Analysis of Per- and Polyfluoroalkyl Substances (PFAS) in Accordance with EPA 1633 Part 2: Analysis of Aqueous Matrices. Waters Application Note [720008143](#). 2023 December.
9. Organtini KL, Rosnack KJ, Hancock P. Expanding the Range of PFAS in a Single Injection to Include Ultra Short Chains Using the Atlantis BEH C18AX Mixed Mode Column. Waters Application Note [720008034](#). 2023 August.
10. Twohig M, Fujimoto G, Mohan A, Organtini K, Rosnack KJ, Hird S. Approaches to Non-Targeted Analyses of Per- and Polyfluoroalkyl Substances (PFAS) in Environmental Samples. Waters Application Note [720007184](#). 2021 March.

11. Charbonnet JA, McDonough CA, Xiao F, Schwichtenberg T, Cao D, Kaserzon S, Thomas KV, Dewapriya P, Place BJ, Schymanski EL, Field JA, Helbling DE, Higgins, CP. Communicating Confidence of Per- and Polyfluoroalkyl Substance Identification via High-Resolution Mass Spectrometry. *Environ.Sci.Technol.Lett* .2022, 9, 473–481
12. Mullin L, Jobst K, DiLorenzo RA, Plumb R, Reiner EJ, Yeung LWY, Jogsten IE. Liquid chromatography-ion mobility-high resolution mass spectrometry for analysis of pollutants in indoor dust: Identification and predictive capabilities. *Analytica Chimica Acta*. 2020 Aug 1125: 29–40
13. Foster M, Rainey M, Watson C, Dodds JN, Kirkwood KI, Fernandez FM, Baker ES. Uncovering PFAS and Other Xenobiotics in the Dark Metabolome Using Ion Mobility, Mass Defect Analysis, and Machine Learning. *Environ.Sci.Technol*. 2022, 56, 12, 9133–9143
14. MacNeil A, Li X, Amiri R, Muir DCG, Simpson A, Simpson MJ, Dorman FL, Jobst KJ. Gas Chromatography- (Cyclic) Ion Mobility Mass Spectrometry: A Novel Platform for the Discovery of Unknown Per-/Polyfluoroalkyl Substances. *Anal.Chem*. 2022, 94: 11096–11103
15. Huset CA, Barlaz MA, Barofsky DF, Field JA. Quantitative Determination of Fluorochemicals in Municipal Landfill Leachates. *Chemosphere*. 2011 Mar 82, 10, 1380–1386
16. Newton S, McMahan R, Stoeckel JA, Chislock M, Lindstrom A, Strynar M. Novel Polyfluorinated Compounds Identified Using High Resolution Mass Spectrometry Downstream of Manufacturing Facilities near Decatur, Alabama. *Environ.Sci.Technol*. 2017, 51, 1544–1552

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© 2025 Waters Corporation. All Rights Reserved.

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